

Release Notes

# GrainMapper3D™ 3.0

Non-destructive 3D Grain Mapping Solution for  
Laboratory Diffraction Contrast Tomography



(This Page Intentionally Left Blank)

# Contents

<b>New Features</b>	<b>4</b>
Projection Geometry Support . . . . .	4
Advanced Acquisition . . . . .	5
DCT Acquisition Wizard . . . . .	6
DCT Data Import with Preview . . . . .	9
DCT Signal and Illumination Regions . . . . .	12
Absorption Data Import with Preview . . . . .	15
<b>Product Enhancements</b>	<b>17</b>
Preview of Absorption Segmentation . . . . .	17
Other Enhancements . . . . .	17
<b>User Interface Changes</b>	<b>18</b>


# New Features

## Projection Geometry Support

GrainMapper3D 3.0 introduces the reconstruction of Flat Panel data collected in projection geometry on the new crystallographic micro-computed tomography system ZEISS Xradia CrystalCT.

Figure 1 displays the geometry setup with example projections for the two supported data collection modes in Laue focusing and projection geometry using the corresponding detectors DCT4X and Flat Panel, respectively. Typical working distances are  $12\text{ mm} < L_{SS} = L_{SD} < 25\text{ mm}$  for Laue focusing geometry and  $10\text{ mm} < L_{SS} < 100\text{ mm}$ ,  $100\text{ mm} < L_{SD} < 550\text{ mm}$  for projection geometry.


**Figure 1** LabDCT Setup for Laue Focusing and Projection Geometries

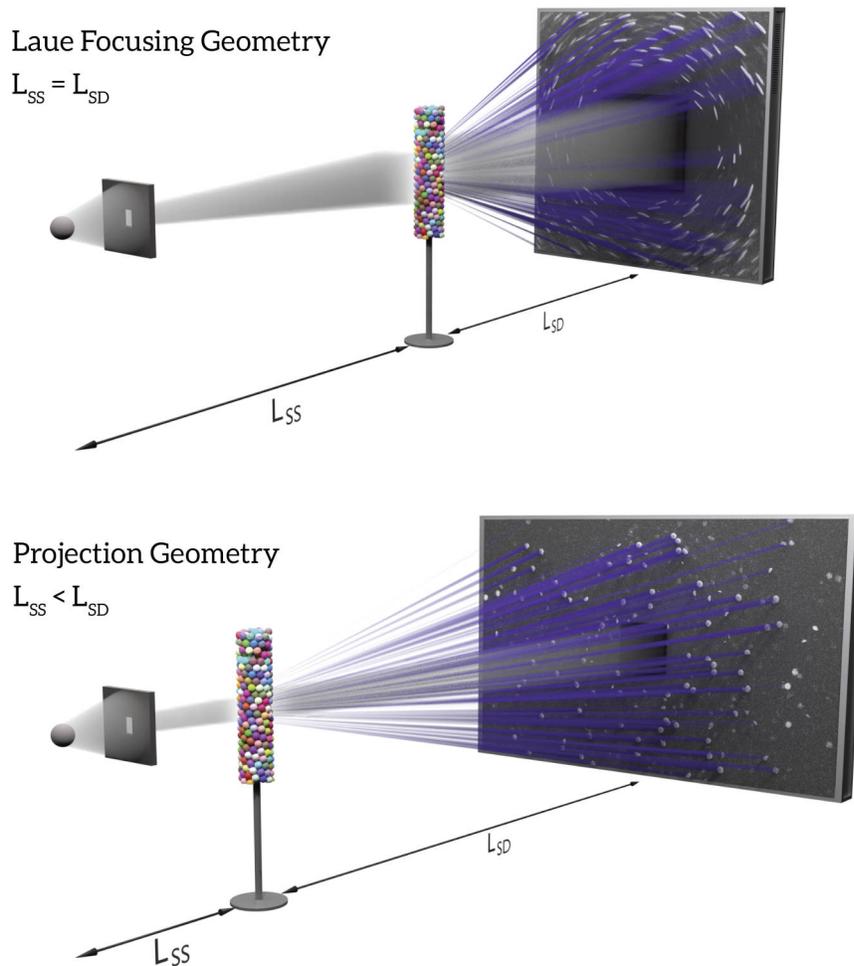

## Advanced Acquisition

Advanced Acquisition allows for seamless collection and subsequent reconstruction of DCT data for larger, irregularly shaped sample volumes as illustrated artistically in Figure 2. Setting up an advanced acquisition of DCT data is done using the new, dedicated DCT Acquisition Wizard, which is described in the next section.

**Figure 2** Advanced Acquisition for Large, Irregularly Shaped Specimens

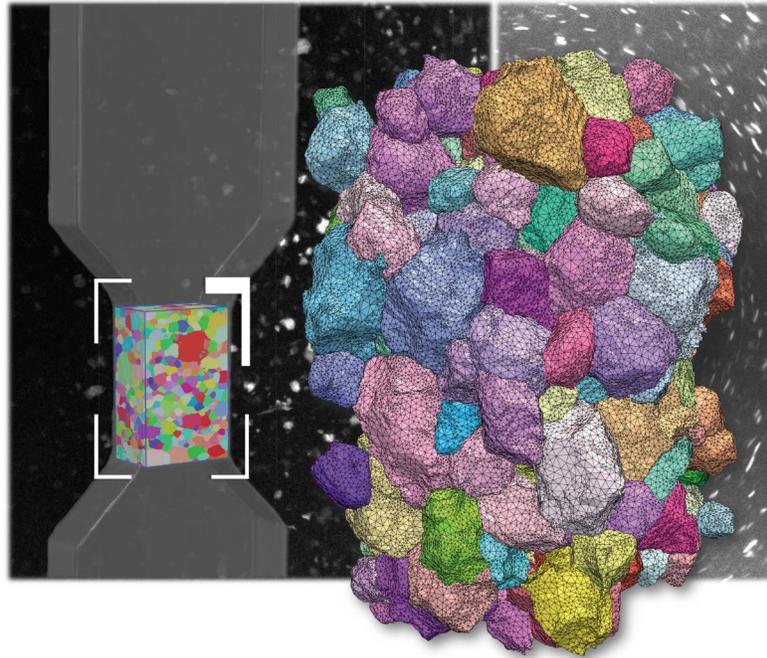

While the *conventional* DCT scan performed using Scout-and-Scan is applicable when the *sample region of interest* (ROI) fits in the *beam field of view* (FOV), the DCT Acquisition Wizard allows for collecting DCT data with three new *advanced* acquisition strategies for samples where the ROI does not fit in the FOV.


Depending on how the ROI exceeds the FOV one can choose between:

- *Helical Phyllotaxis* covers samples with a vertical extent larger than the beam height by means of a “Golden Angle” rotation of  $\sim 137.5^\circ$  combined with a vertical translation of the order  $\sim 1\text{--}5\ \mu\text{m}$  between consecutive projections.
- *Helical Phyllotaxis Raster* is used if the ROI exceeds the FOV both vertically and horizontally. It combines the helical phyllotaxis rotation and vertical translation with a horizontal translation. The number of horizontal steps is the same for all projection angles.
- *Helical Phyllotaxis HART* (high aspect ratio tomography) is tailored for specimens with plate-like geometries. It combines the helical phyllotaxis rotation and vertical translation with a horizontal translation. The number of horizontal steps is adapted to tightly fit the ROI at every projection angle.

The advanced acquisition strategies are schematically compared to the conventional DCT scan in Figure 3. Here the points, colored by the sample rotation angle, mark where the center of the beam intersects the sample surface for individual projections.

**Figure 3** Advanced Acquisition Strategies for Different Specimen Shapes

Scanning modes	Conventional	Helical Phyllotaxis	Helical Phyllotaxis Raster	Helical Phyllotaxis HART
FOV vs. ROI	ROI fits in FOV	ROI taller than FOV	ROI larger than FOV	ROI larger than FOV
Rotation stepping	$360^\circ/N$	$137.5^\circ$	$137.5^\circ$	$137.5^\circ$
Vertical translation	NO	YES	YES	YES
Horizontal translation	NO	NO	YES	YES - adaptive

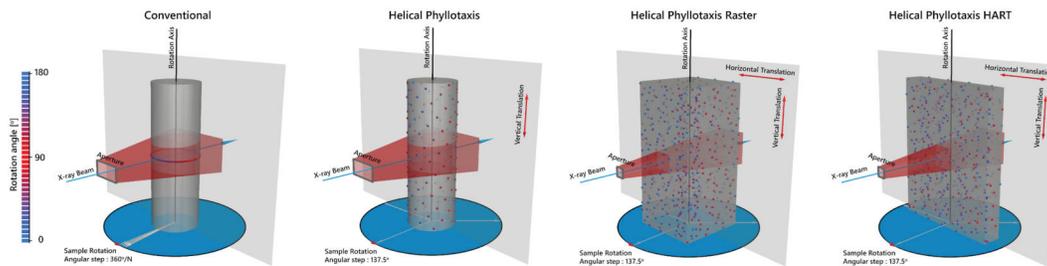

## DCT Acquisition Wizard

The **DCT Acquisition Wizard** is a new tool to help set up an advanced acquisition DCT scan. Only the principal structure and a few selected features of the DCT Acquisition Wizard will be outlined here.



The reader is referred to the GrainMapper3D User's Guide 3.0 for full instructions.




The DCT Acquisition Wizard is only to be installed on the acquisition workstation by ZEISS Service.

In order to perform an advanced acquisition of DCT data with the DCT Acquisition Wizard, select **Start > All Programs > Xnovo Technology ApS > Xnovo Acquisition Wizard** from the **Windows Start** menu. The **DCT Acquisition Wizard** window opens and requests the exclusive control over the Scout-and-Scan control system.

Figure 4 shows the layout of the DCT Acquisition Wizard, organized in **Sample** section, **Acquisition** section (including tabs for **Acquisition Settings**, **Acquisition Strategy**, **Acquisition Options**, and **Sample Position**), and **Recipe** section.


**Figure 4** Principal Layout of the DCT Acquisition Wizard

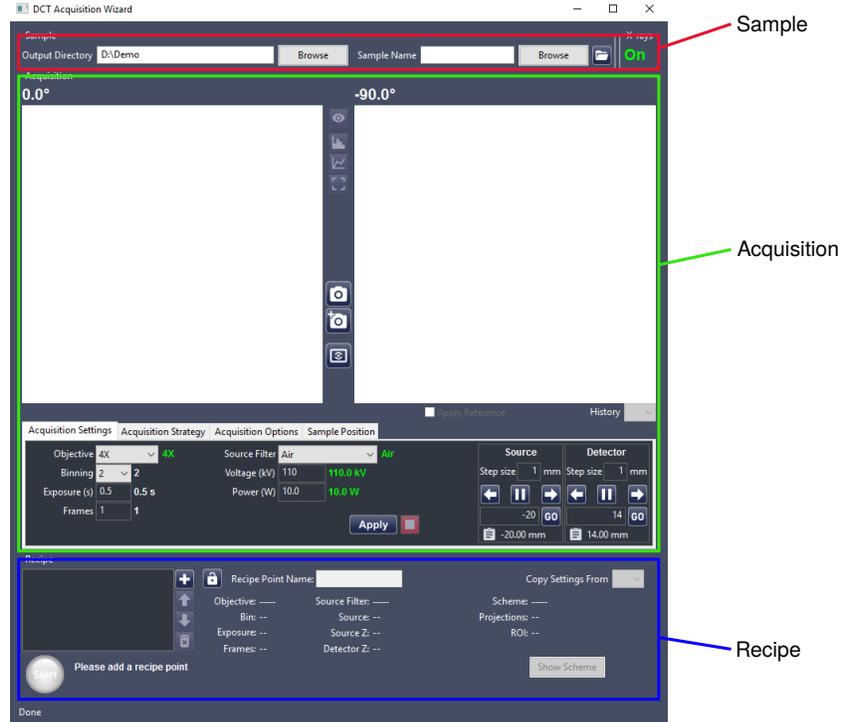

The *look and feel* of the DCT Acquisition Wizard is similar to Scout-and-Scan, and the user should be able to recognize many of the controls, for instance in the **Acquisition Settings**, **Sample Position** and **Recipe** sections.


**Defining the ROI** is straightforward in the DCT Acquisition Wizard. Irrespective of the beam FOV, simply drag rectangles on top of the  $0^\circ$  and  $-90^\circ$  absorption projections to cover the desired sample ROI as illustrated in [Figure 5](#). The figure also shows the *FOV coverage* with respect to the ROI as the partly overlapping red boxes corresponding to the selected acquisition settings and strategy.

**Figure 5** Defining the ROI in the DCT Acquisition Wizard

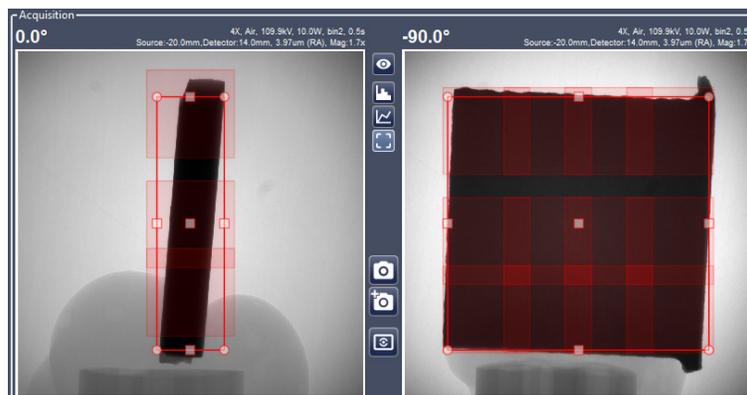

The new *advanced* acquisition strategies outlined in [Figure 3](#) can be selected in the **Strategy** drop-down of the Acquisition Strategy tab, see [Figure 6](#).


**Figure 6** Defining the Acquisition Strategy in the DCT Acquisition Wizard

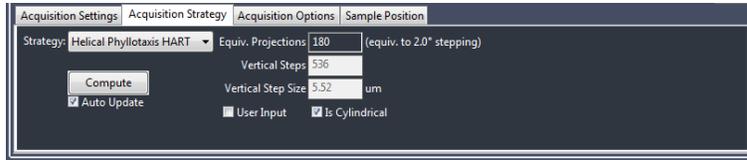

The number of **Equivalent Projections** must be provided by the user based on the detector working distance. The scanning modes with Helical Phyllotaxis motion then fits this number of projections into the aperture FOV height by adjusting the **Vertical Step Size** between consecutive images accordingly. The number of **Vertical Steps** then follows to cover the entire height of the ROI defined by the user. Uncheck **Is Cylindrical** if applicable for the selected acquisition strategy and the sample has square or rectangular shape.

[Figure 7](#) show the DCT Acquisition Wizard during a typical scan. The progressing scan views the latest projection in the Acquisition section, while the estimated time remaining is given next to the progress bar at the bottom.


**Figure 7** DCT Scan in Progress in the DCT Acquisition Wizard

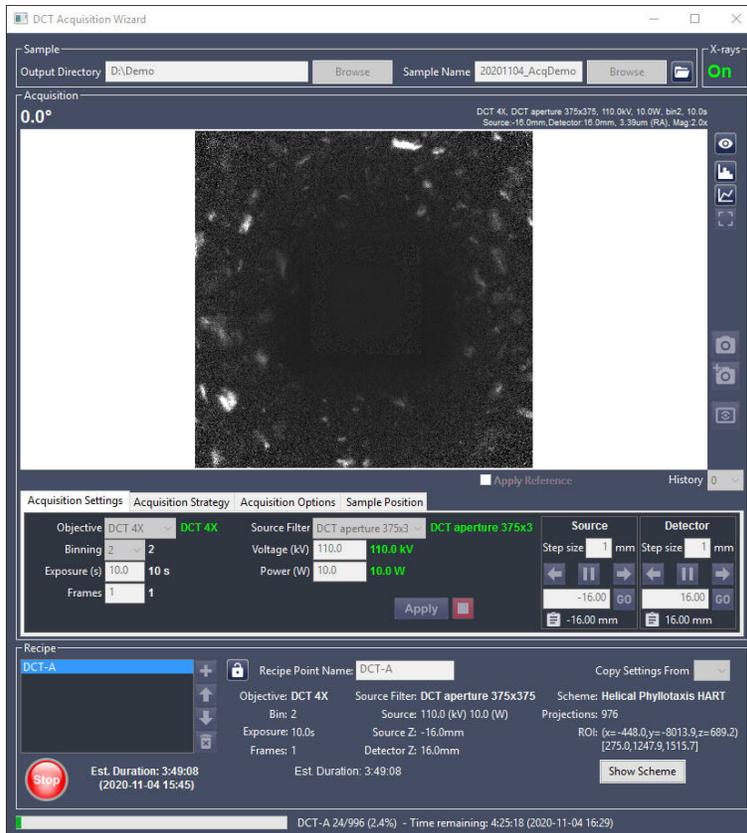

## DCT Data Import with Preview

The **Project** tab of GrainMapper3D has been redesigned to make space for a more sophisticated **DCT Data** import section with a live **DCT Data Preview**. The **Project** tab now focuses on the choice of appropriate DCT data import parameters to perform a better background correction with feedback. The absorption data import has been moved to the **Absorption Mask** tab. [Figure 8](#) shows the redesigned **Project** tab for a new project organized in sections **Sample**, **DCT Data** and **Layers**. Parts of the former advanced DCT import options have been migrated into a new **Background Correction** group and **Noise Reduction** group.

**Figure 8** New Layout of Project Tab with DCT Data Import Preview

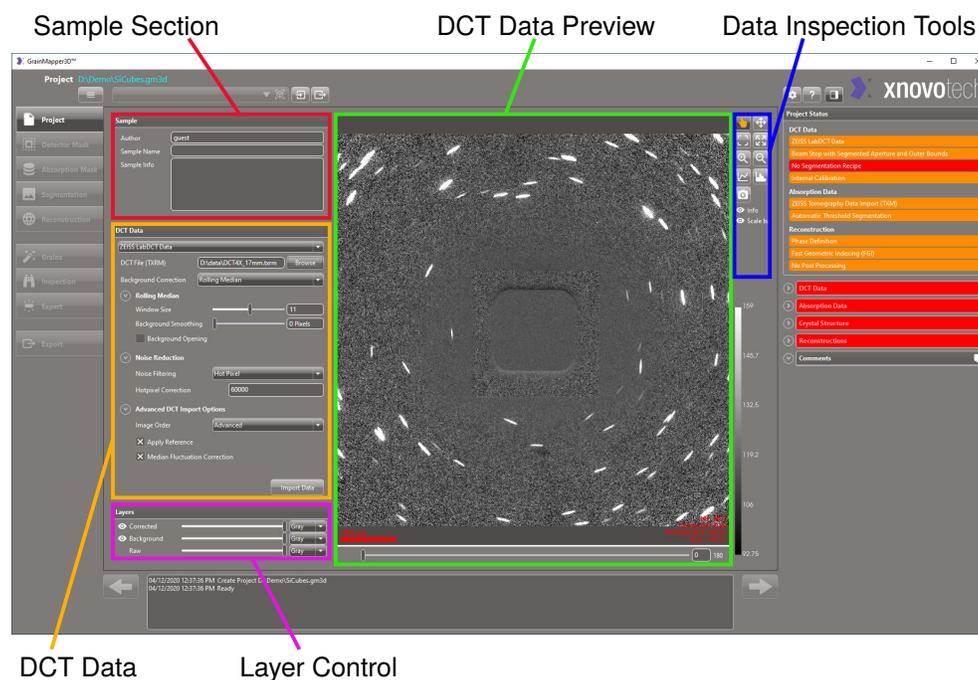

**Background Correction** The former DCT data import options have been moved into the **Rolling Median** background correction method and a new **Iterative Background** correction method has been added. [Table 1](#) details the new controls for **Background Correction** of the Diffraction Contrast Patterns (DCPs).



**Noise Reduction** A new noise reduction method **Denoising** has been added. [Table 2](#) gives an overview of the new controls for **Noise Reduction** of the DCPs. The noise reduction is performed on the raw DCPs after reference correction and prior to background calculation and subtraction.

**Advanced Acquisition Support** By default, the DCPs are resorted such that similar sample positions follow each other. Sorting is important for samples with a footprint larger than the beam stop, where sample attenuation contributes to the background pattern depending on both scan angle and sample position. Sometimes, mostly for


helical phyllotaxis scans without horizontal sample translations, there are artifacts in the background that arise from the sample absorption and can be reduced by changing the **Image Order** to **Scan Order**, see [Table 3](#).

**Layer Control** The visual appearance of the preview can be adjusted using the **Layer** control as described in [Table 4](#). A cache for the DCT Data Import Preview allows quicker switching back and forth between already calculated import settings.

**Table 1** New Controls for DCT Background Correction in Project Tab

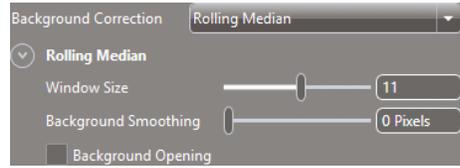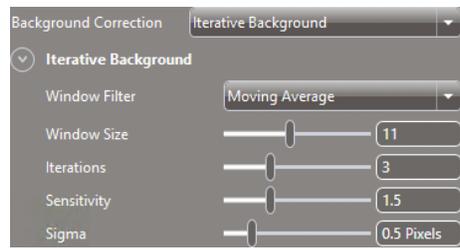Control	Function
	<p><b>Rolling Median Background Correction for DCT Data:</b></p> <ul style="list-style-type: none"> <li>• <b>Window Size</b>, apply the rolling median correction over the specified number of successive images. Subtracts pixel-wise median of image stack.</li> <li>• <b>Background Smoothing</b>, apply a 2D Gaussian filter to the rolling median background image with the specified sigma value (in pixels). Should only be used if no reference was collected.</li> <li>• <b>Background Opening</b>, check to apply a morphological opening denoising step on the rolling median background image.</li> </ul>
	<p><b>Iterative Background Correction for DCT Data:</b></p> <ul style="list-style-type: none"> <li>• <b>Window Filter</b>, choose between: <ul style="list-style-type: none"> <li>– <b>Moving Average</b>, average over images in window,</li> <li>– <b>Rolling Median</b>, take median over images in window, or</li> <li>– <b>Gaussian</b>, images in window will be averaged with Gaussian weighting.</li> </ul> </li> <li>• <b>Window Size</b>, apply the above selected window filter over the specified number of images.</li> <li>• <b>Iterations</b>, specify the number of adaptive background interpolation steps to be performed.</li> <li>• <b>Sensitivity</b>, set intensity threshold over window average, above which intensity is replaced by interpolated value.</li> <li>• <b>Sigma</b>, set Gaussian shape parameter for 2D smoothing at every iteration.</li> </ul>

**Table 2** New Controls for DCT Noise Reduction in Project Tab

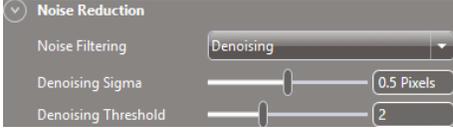Control	Function
	<p><b>Noise Reduction for DCT Data</b>, choose between the following <b>Noise Filtering</b> options:</p> <ul style="list-style-type: none"> <li>• <b>Hot Pixel</b>, set maximum allowed intensity. Higher values are assumed to be outliers and sets to zero,</li> </ul>
	<ul style="list-style-type: none"> <li>• <b>Denoising</b> according to: <ul style="list-style-type: none"> <li>– <b>Denoising Sigma</b>, set 2D Gaussian shape parameter for calculating the smoothed background noise level.</li> <li>– <b>Denoising Threshold</b>, set threshold over background noise level, above which intensity is replaced by interpolated value.</li> </ul> </li> </ul>
	<ul style="list-style-type: none"> <li>• <b>None</b>, to not apply any noise filtering.</li> </ul>

**Table 3** New Control for the Image Order in Project Tab

Control	Function
	<p><b>Advanced DCT Import Options for DCT Data:</b></p> <ul style="list-style-type: none"> <li>• Change <b>Image Order</b> to <b>Scan Order</b> if there are sample absorption artifacts in the images that are better reduced by this image order. Usually only applicable for helical phyllotaxis scans.</li> </ul>

**Table 4** Layer Control for the DCT Data Import Preview in Project Tab


Control	Function
	<p><b>Layers</b> is the control for visualizing the different corrections applied to the imported <b>DCT Data</b>:</p> <ul style="list-style-type: none"> <li>• <b>Raw</b> DCPs (bottom, always shown)</li> <li>• <b>Background</b> images for each Raw DCP (middle),</li> <li>• <b>Corrected</b> DCPs (top)</li> </ul> <p>Use the eye to turn visibility on/off.  Use the sliders to change opacity of the layers.  Use the drop down boxes to change colors of the layers.</p>

## DCT Signal and Illumination Regions

The **Detector Mask** tab is now the second processing step of the GrainMapper3D workflow. The order has been changed in order to exploit the new *aperture illumination* feature when defining the ROI during the absorption data import in the subsequent **Absorption Mask** tab.


Figure 9 shows the revised layout of the **Detector Mask** tab organized into a data processing **Detector Mask** section, **Layers** control, **DCT Data View** with **Data Inspection Tools**.

**Figure 9** DCT Signal and Illumination Regions in Detector Mask Tab

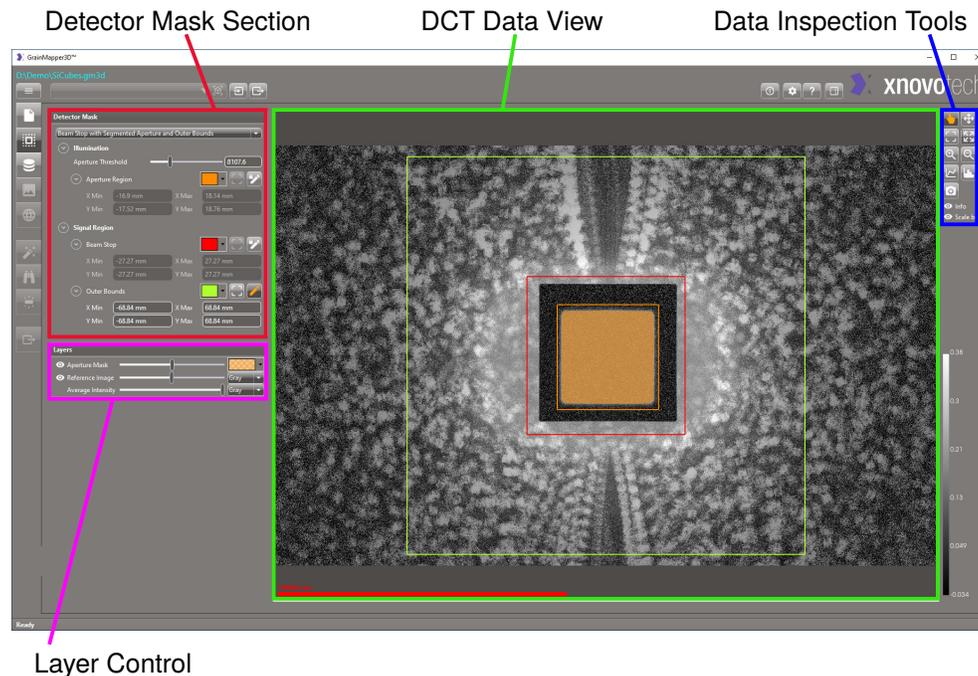





The **DCT Data View** now also shows the **Reference Image** if present. The visual appearance can be adjusted using the **Layers** control as described in Table 5.

**Table 5** Layer Controls of the Detector Mask Tab

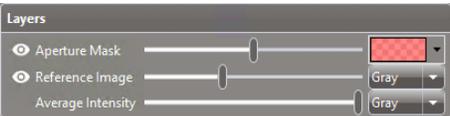Control	Function
	<p>Use the <b>Layers</b> control to visualize:</p> <ul style="list-style-type: none"> <li>• <b>Average Intensity</b> of all DCPs (bottom, always shown)</li> <li>• <b>Reference Image</b> (middle)</li> <li>• <b>Aperture Mask</b> (top – applies only to <b>Segmented Aperture</b>)</li> </ul> <p>Use the eye to turn visibility on/off.            Use the sliders to change opacity of the layers.            Use the drop down boxes to change colors of the layers.</p>

The first step on the **Detector Mask** tab is to choose the appropriate **Detector Mask Recipe**. The list of detector mask recipes has been revised and the recipes with predefined beam stop sizes have been replaced. The available recipes are listed in [Table 6](#) along with a description of when to use which recipe. Start from the top of the table and choose the first recipe that fits the current DCT Data.

**Table 6** Controls of Detector Mask Tab

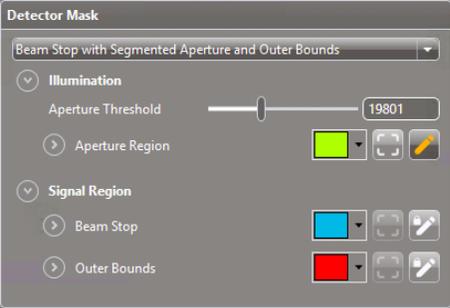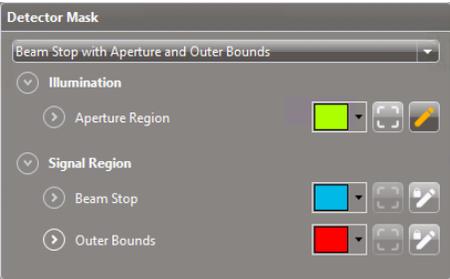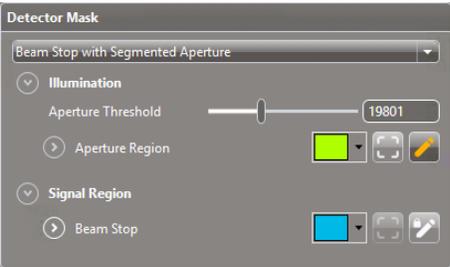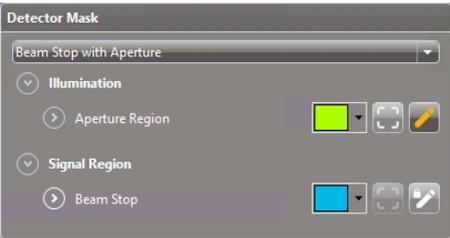Control	Function
	<p><b>Beam Stop with Segmented Aperture and Outer Bounds</b>, use if:</p> <ul style="list-style-type: none"> <li>• The aperture illumination is visible through the area covered by the beam stop</li> <li>• The detector signal towards the edges is noisy and/or very weak</li> </ul>
	<p><b>Beam Stop with Aperture and Outer Bounds</b>, use if:</p> <ul style="list-style-type: none"> <li>• The automatic aperture segmentation is not applicable</li> <li>• The detector signal towards the edges is noisy and/or very weak</li> </ul>
	<p><b>Beam Stop with Segmented Aperture</b>, use if:</p> <ul style="list-style-type: none"> <li>• The aperture illumination is visible through the area covered by the beam stop</li> <li>• The detector signal is good all the way to the edges</li> </ul>
	<p><b>Beam Stop with Aperture</b>, use if:</p> <ul style="list-style-type: none"> <li>• The automatic aperture segmentation is not applicable</li> <li>• The detector signal is good all the way to the edges</li> </ul>
	<p><b>Beam Stop</b>, use only for:</p> <ul style="list-style-type: none"> <li>• <i>Conventional</i> DCT data where it is not necessary to define the aperture illumination</li> </ul>


**Illumination – Aperture Region** For any DCT scan where the sample is larger than the aperture FOV, a recipe allowing the user to set the region of direct illumination, defined by the aperture, must be selected. This is done in order to correctly compute which part of the sample is illuminated.

For any *advanced* DCT scan choose a recipe defining the aperture region. It is also recommended to use a recipe with aperture illumination definition for *conventional* DCT scans.

The controls to define the aperture can be seen in [Table 7](#). There are two different recipe types for this:

- **Aperture**, which defines the aperture as a centered box of preset size (375x375um, 250x750um or 750x750um) that can be adjusted by dragging or typing in the boundaries
- **Segmented Aperture**, which allows the user to segment the aperture inside the defined box by setting a threshold.

**Table 7** Controls of the Aperture

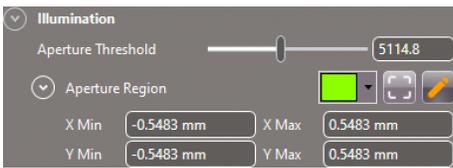Control	Function
	<p>If the sample ROI is larger than the aperture FOV it is important to define the <b>Aperture Illumination</b>, either:</p> <ul style="list-style-type: none"> <li>• Drag the <b>Aperture Threshold</b> slider to set the segmentation threshold for recipes with <b>Segmented Aperture</b></li> <li>• Press  to select one of the predefined apertures</li> <li>• Press  to edit the aperture FOV</li> <li>• Give the boundaries of the aperture</li> </ul>

**Signal Region – Beam Stop** The region on the detector, which is not hit by diffraction spots due to the presence of the beam stop, must be masked out. This must be done for all detector mask recipes.

The beam stop used during acquisition needs to be entered. It is not carried over in the DCT file metadata, as it is mounted manually onto the detector. The predefined beam stops will appear centered, but can be moved and adjusted in size.

The controls to define the beam stop can be seen in [Table 8](#).

**Signal Region – Outer Bounds** Using a detector mask recipe with outer bounds is recommended if the detector signal towards the edges is weak or noisy, for instance due to self-absorption. This usually applies to DCT data collected with a Flat Panel detector.

**Table 8** Controls of the Beam Stop

Control	Function
	<p>Define the <b>Signal Region</b> covered by the <b>Beam Stop</b>, either:</p> <ul style="list-style-type: none"> <li>• Press  to select one of the preset beam stop sizes: <ul style="list-style-type: none"> <li>– For the DCT 4X objective</li> <li>– For the Flat Panel detector</li> </ul> </li> <li>• Press  to edit the beam stop</li> <li>• Give the boundaries of the beam stop</li> </ul>

**Table 9** Controls of the Outer Bounds

Control	Function
	<p>Define the <b>Outer Bounds</b> of the <b>Signal Region</b> if the detector signal towards the edges is noisy and/or very weak, either:</p> <ul style="list-style-type: none"> <li>• Press  for preset values for the Flat Panel detector: <ul style="list-style-type: none"> <li>– Default Outer Bound</li> <li>– Outer Bounds enclosing 50, 60, 70, 80, 90, or 100% of the detected intensity</li> </ul> </li> <li>• Press  to edit the outer bounds</li> <li>• Give the boundaries of the outer bounds</li> </ul>

The controls to define the outer bounds can be seen in [Table 9](#). The preset percentile outer bounds are centered on the detector and enclose the named fractions of the detected intensity.


## Absorption Data Import with Preview

The absorption data import, which was previously done in the **Project** tab, has been moved to a dedicated section in the **Absorption Mask** tab. The absorption crop is now done when importing the data. This processing order has the advantage that the *aperture illumination*, defined in the **Detector Mask** tab, now falls prior to the absorption data import, such that the actual illuminated region can be exploited when defining the ROI for the absorption data.

[Figure 10](#) shows the principal layout of the *data import view* of the **Absorption Mask** tab, organized into an **Absorption Data** section for importing data, an **Absorption Data Preview** view with **Data Inspection Tools**, a **Preview** layer control section and a **Vertical ROI Control**.

[Table 10](#) details the control to import absorption data. Once an **Absorption File**

**Figure 10** Absorption Data Import in Absorption Mask Tab

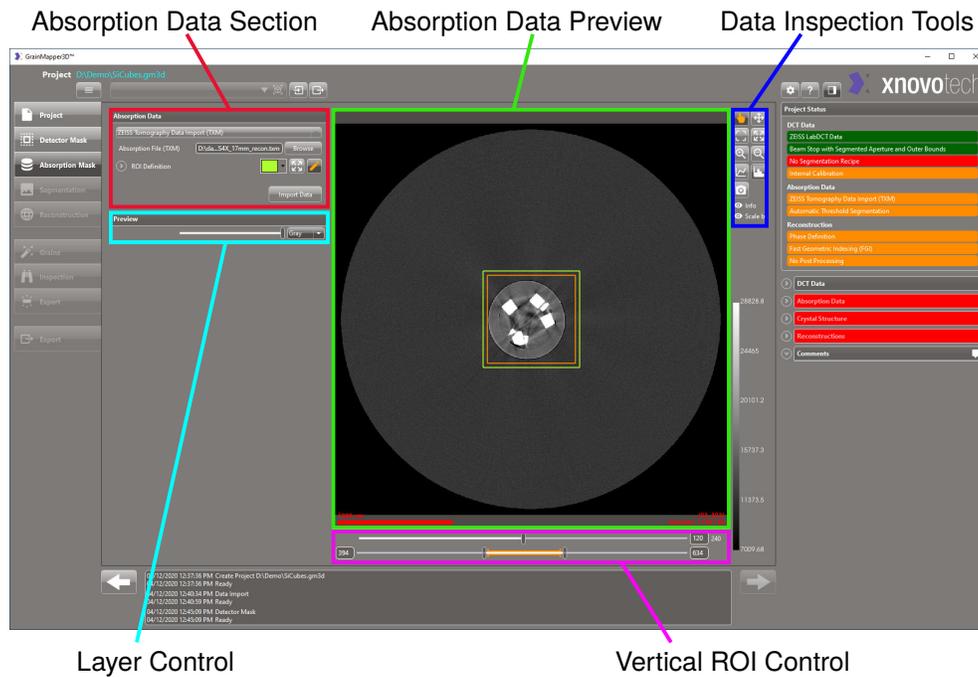


(TXM) has been specified, the **Absorption Data Preview** displays the horizontal slice of the absorption data centered in the aperture FOV. The **orange** rectangle on the preview indicates the horizontal aperture FOV, while the ROI Definition is shown in **green**.

The bottom line of the **Vertical ROI Control** corresponds to all vertical slices in the absorption data, the **orange** line segment corresponds to the vertical aperture FOV and the two sliders correspond to the top and bottom ROI limits, which can also be input as slice numbers. The top slider controls the slice preview between the first and last image of the selected ROI.

Figure 11 shows the principal layout of the **Absorption Mask** tab after importing the absorption data.

**Table 10** Controls of Absorption Mask Tab

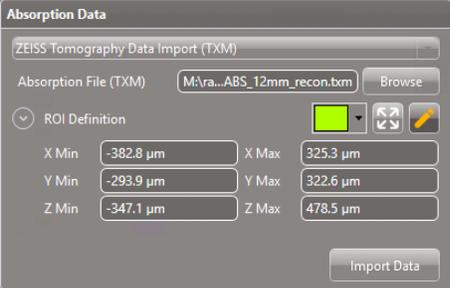
Control	Function
	<p>ZEISS Tomography Data Import (TXM) imports an absorption volume.</p> <p>To locate the Absorption File (TXM), press Browse.</p> <p>ROI Definition specifies the spatial coordinates of the ROI prior to data import.</p> <p>Press  to freely select the ROI (<b>green</b>) relative to the aperture FOV (<b>orange</b>).</p> <p>Press Import Data to import the cropped <i>absorption data</i>.</p>

## Product Enhancements

### Preview of Absorption Segmentation

The absorption **Threshold Segmentation** and **Range Segmentation** recipe now supports live preview of the chosen threshold or range, respectively. Whenever the threshold or range is changed, a preview layer is updated to indicate which values will be masked. The **Preview** layer can be toggled on or off in the Layer control. However, the live preview only segments individual slices and does not support segmentation filters.

**Figure 11** Absorption Data Segmentation in Absorption Mask Tab

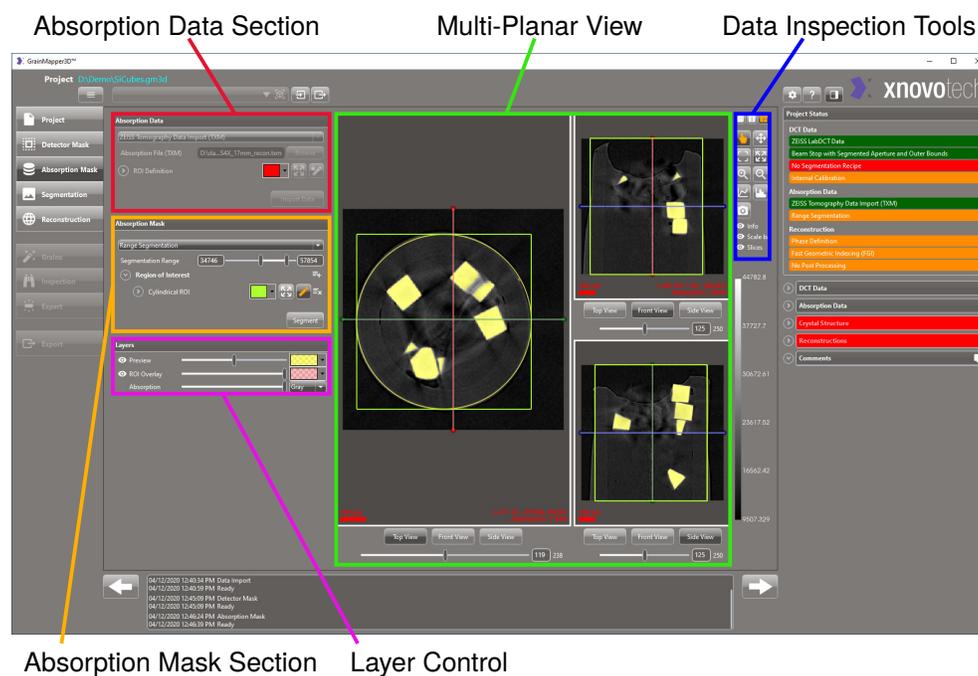

### Other Enhancements

**Importing Combined Segmentation** When importing recipes from another project file, a combined segmentation is now automatically imported together with the depended segmentations without actively selecting them.

**Result File** If the user desires to **Export IPF Color Coding** in the Export tab, it is now possible to choose between two different **Data Types**, namely FLOAT and UINT8. The UINT8 option was added as it corresponds a significant reduction in output file size.

**Visualization** The underlying graphics library has been updated which results in minor rendering improvements.

## User Interface Changes

**Flat Panel Support** Flat Panel data is fully supported in GrainMapper3D 3.0, both as DCT and absorption data. The program automatically detects the detector used and adjusts relevant recipes accordingly, for instance:


- Preset **Aperture**, **Beam Stop** and **Outer Bound** in the **Detector Mask** tab
- **Self Fitting** parameters and previews in the **Expert** tab

**Project File** The project file format has updated in order to adapt to the new features introduced in GrainMapper3D 3.0. This means that a GrainMapper3D 3.0 project file cannot be opened with any prior version of GrainMapper3D. Vice versa, if the user attempts to open a prior GrainMapper3D 2.x project file in GrainMapper3D 3.0 a warning will be issued that the file format is about to change. It is recommended to open prior GrainMapper3D project files as read-only in GrainMapper3D 3.0 to preserve the prior file format or to make a copy\_v3.gm3d before opening.

**Project Status** The recipes listed in the **Project Status** section have been re-arranged into the groups **DCT Data**, **Absorption Data** and **Reconstruction** to accommodate the new processing workflow.

**Histogram control** The histogram controls when importing DCT or absorption data now have a **Global Range** checkbox as shown in [Figure 12](#). Check to preserve the data range when changing processing parameters or slicing through the data sets.

**Figure 12** Histogram Control with Global Range

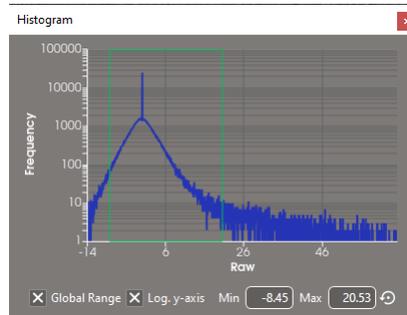

**New Preferences** A few additional preferences have been added, allowing the user to tailor the new recipes in the user interface, namely:

- Preferred **Background Correction** and **Noise Reduction** recipes for **DCT Data Import**.
- Preferred **Data Type** for **Export IPF Color Coding**.

(This Page Intentionally Left Blank)

## **GrainMapper3D™**

Release Notes

### **Xnovo Technology ApS**

Galoche Alle 15, 1.

4600 Køge, Denmark (DK)

support@xnovotech.com

www.xnovotech.com

EAN: DK-34894221

© 2021 by Xnovo Technology ApS

This document or any part of it must not be translated, reproduced, or transmitted in any form or by any means, electronic or mechanical, including photocopying, recording, or by any information or retrieval system. Violations will be prosecuted.

The use of general descriptive names, registered names, trademarks, etc. in this document does not imply, even in the absence of a specific statement, that such names are exempt from the relevant protective laws and regulations and therefore free for general use. Software programs will fully remain the property of Xnovo Technology ApS. No program, documentation, or subsequent upgrade thereof may be disclosed to any third party, unless prior written consent of Xnovo Technology ApS has been procured to do so, nor may be copied or otherwise duplicated, even for the customer's internal needs apart from a single back-up copy for safety purposes.

Xnovo Technology ApS reserves the right to make modifications to this document without notice.